


ATUMtome Pre-installation Considerations

RMC
Products
by Boeckeler®

BOECKELER INSTRUMENTS, INC.
4650 S. Butterfield Drive
Tucson, Arizona 85714 USA
Phone: 520.745.0001
Fax: 520.745.0004
rmcproducts.com

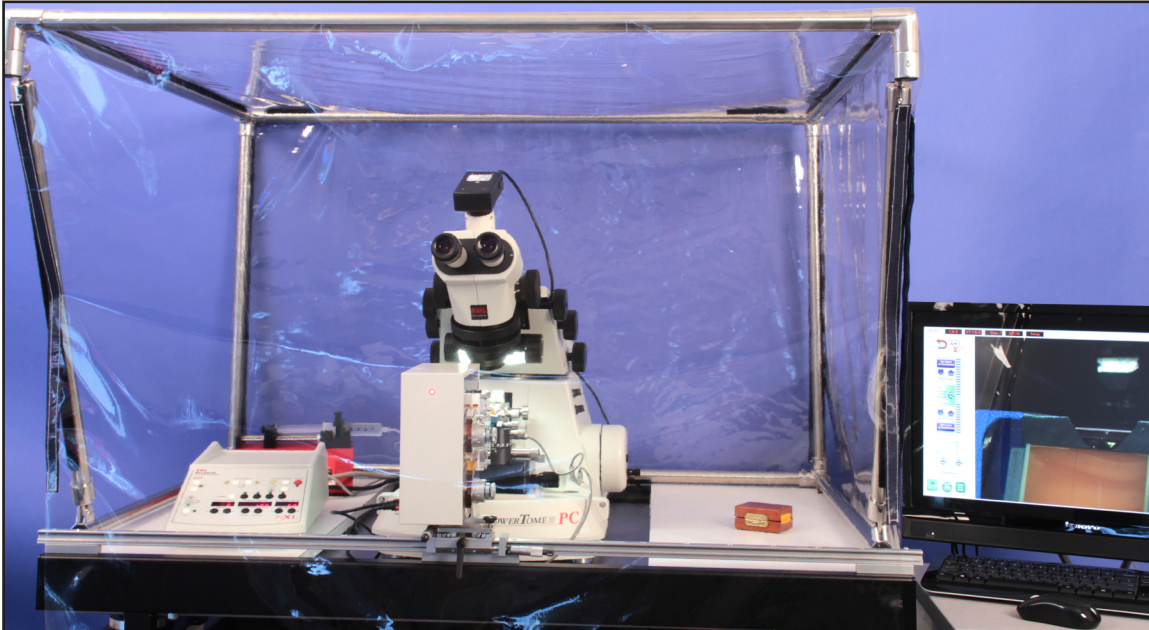

Location (room): suggest 8ft x 8ft (2.5m x 2.5m) minimum
Free of air drafts (draughts) and temperature fluctuations. Air conditioning ducts should be located to avoid directing air onto the equipment.

Humidity: Relative humidity ideally 45%rh

Doorway opening: minimum 32in (81cm)

Floor must bear weight of table and ultramicrotome: 839lbs (381kg)

Space for air table: 48in x 30in (122cm x 76cm)

Space for computer: minimum 26in x 24in (66cm x 61cm)

Water required for pump: double distilled (30ml per run)

Air or Nitrogen supply: 90-100PSI (6-7bar)

Electricity: 120V 10A or 240V 5A (located within 3ft (1m))

Internet: needed for tech support through Skype